

GermBank

Bacterial strain preservation system

Features:

- Easy to use system simplify the storage and retrieval steps of bacterial cultures
- Each vial contains 25 beads with cryopreservative ensures longer survival of fastidious bacteria
- 5 colors of tubes are easy distinction



Preservation steps

1. inoculate colonies from a fresh, pure culture.
2. Cap back tightly and invert the tube 4-5 times.
3. Remove excess cryopreservative using sterile pipette.
4. Store in appropriate freezer or liquid nitrogen.

Recovery steps

1. Insert a sterile needle through the hole.
2. The bead may then be used to inoculate appropriate solid or liquid media.

